

SanAir Dust & Mite Collector



- Collection and Extraction within one single device
- No dust handling necessary: Only 2 minutes sampling time
- One Collector per Sample
- Allows exposure to be expressed per weight or per unit area
- Cost effective

The Mitest collector is a moulded plastic device with an inner dust/mite collector which fits onto the wand of a vacuum cleaner. After collection, a cap is applied to the bottom of the collector.

Dust samples can be collected from bedding, soft furnishings or carpets and extracted within 10 minutes.

Dust sampling using the Mitest Collector

Sampling Site	n	mg of dust collected (mean +/- SD)
Bed/Mattress	12	140 +/- 85
Carpet	27	660 +/- 446
Sofa	7	142 +/- 151

Instructions for Use of the Mitest Collector

Attach the collector to the vacuum cleaner tube. Turn on the vacuum cleaner and vacuum 4 separate 8.5 x 11" areas (letter-size paper) for 30 seconds for each suspect item (chair, carpet etc.) Use one collector cassette for each sample area. Use one collector/cassette in a non-suspect area as a control sample.